EN	ES	IT	NL
DA	NO	AR	JP
FR	DE	PT	EL
FI	SV	RU	UA
CS	HU	LV	PL
BG	SL	KOR	TH
ET	LT	MT	RO
SK	TR	MS	ZH

Single Use Garment Type 5/6

INSTRUCTIONS FOR USE

Ansell

Single Use Garment Type 5/6

This Instruction for Use note is to be used in combination with the specific information that is mentioned on each packaging enclosure.

USE OF GARMENT:

This is a single use cleanroom garment designed to provide product protection and which protects against fine dry particles (Type 5) and limited liquid splashes and fine sprays (Type 6). Check the garment carefully to ensure there are no defects. Ensure the expiry date has not been reached.

For full protection wear compatible gloves, facemask, goggles and footwear.

WEARING GARMENT:

Don garment, zip up fully and press zipper into lock position. Remove backing tape from inside of zip cover and seal zip cover to garment - ensuring zip is fully covered.

For full protection wear compatible gloves, facemask, goggles and footwear offering at least equivalent protection. These should be taped together using waterproof tape to cover each join.

LIMITATIONS OF USE:

Not suitable for prolonged or intensive exposure to chemical spray and splash. Heat stress should not occur but consideration should be given to alternation of work and rest periods. The user is the sole judge of the suitability of the garment, and any accessories, given its intended use and duration.

Nitritex Ltd accepts no responsibility for improper use.

GARMENT FABRIC:

Polyethylene/Polypropylene laminate.

WARNING:

The person wearing the electrostatic dissipative protective clothing shall be properly earthed. The resistance between the person and the earth shall be less than 10⁸ Ω.

Electrostatic dissipative protective clothing shall not be open or removed whilst in presence of flammable or explosive atmospheres or while handling flammable or explosive substances;

Electrostatic dissipative protective clothing shall not be used in oxygen enriched atmospheres without prior approval of the responsible safety engineer; The electrostatic dissipative performance of the electrostatic dissipative protective clothing can be affected by wear and tear, laundering and possible contamination.

Electrostatic dissipative protective clothing shall permanently cover all non-complying materials during normal use (including bending and movements).

STORAGE & TRANSPORT:

Store in a clean, dry place away from heat and light. To prevent the risk of damage, transport in cardboard box provided and retain unopened until required.

DISPOSAL:

Remove carefully and turn inside out to trap contamination. Assess any contamination that may have been introduced during use and destroy immediately in accordance with regulations.

To the best knowledge of the manufacturer, BioClean single use garments do not contain any constituent substances which could lead to skin irritations or skin diseases or other allergic reactions or that are carcinogenic, toxic to reproduction or mutagenous.

CAUTION:


Keep plastic bag away from children to prevent the danger of suffocation.

EXPLANATION OF GARMENT LABEL - KEY:

1. Overall Manufacturer
2. Trademark
3. Model Identification
4. CE marking - complies with Regulation (EU) 2016/425 relating to Category III PPE.
Under the supervision of Notified body - SGS Fimko Oy, P.O. Box 30 (Särkinimestie 3), 00211 HELSINKI, Finland (EU Notified Body No. 0598).
5. Wearer should read these instructions
6. Indicates compliance with EU standards for protective clothing: Type 5 standard EN ISO 13982-1:2004+A1:2010 for fine dry particles; Type 6 standard EN 13034:2005+A1:2009 for reduced spray, limited splash & EN1149-5:2008 antistatic.

7. International Care Symbols a) Do not wash b) Do not chlorine bleach c) Do not tumble dry d) Do not iron e) Do not dry clean f) Inflammable – keep away from naked flame and intensive heat
8. Garment size - Please ensure that you select the correct size of garment according to the SIZING TABLE

9. DO NOT REUSE

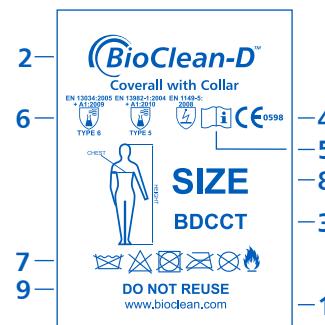The EU Declaration of conformity may be downloaded from:
www.ansell.com/regulatory

EU Type examination carried out by:
Centexbel-Gent, Technologiepark 70, BE-9052 Zwijnaarde (Gent), Belgium.
Notified Body Number 0493

TESTING

The BioClean-D range of PPE Category III Type 5/6 Single use garments have been subject to the following comprehensive tests:

Fabric Property Tests	Test Method	Class or Result
Abrasion Resistance	EN 530 Method 2	1
Trapezoidal Tear Resistance machine dir./x dir.	ISO 9073-4	3/2 (55.5N/29.3N)
Flex Cracking Resistance	ISO 7854 Method B	3
Seam Strength	ISO 13935-2	2 (70N)
Puncture Resistance	EN 863	1 (8N)
Tensile Strength Machine dir./x dir.	EN ISO 13934-1:1999	2/1 (97N/48N)
Charge Decay Time	EN 1149-3: 2004	0.07 sec
Particle Shedding	IEST-RP-CC-003.4 2011 (Helmke Drum)	1

Resistance to Penetration by Liquids EN 6530:2005

Liquid	Repellency Index	Penetration Index
Sulphuric Acid (30%)	3	3
Sodium Hydroxide (10%)	3	3
o-Xylene	2	3
Butan-1-ol	3	3

Whole Suit Test Performance	Test Method	Result
Reduced Spray Test (Type 6)	EN 13034:2005+A1:2009 & EN 468 modified*	Pass
Whole Suit Inward Leakage (Type 5)	EN/ISO 13982-1:2004+A1:2010*	$L_{p90} \text{ 82/90 } < 30\% L_s, 8/10 < 15\%$

*All interfaces with additional PPE were taped during testing, as required by the standards.

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IMPORTER

UK

Nitrilex Ltd
Minton Enterprise Park,
Oaks Drive, Newmarket, Suffolk
CB8 7YY, UK

T: +44 1638 663338
F: +44 1638 668890

EU

Ansell Healthcare Europe NV,
Riverside Business Park,
Blvd International, 55
1070 Brussels, Belgium

T: +32 2 528 7400
F: +32 2 528 7401

Australia

Ansell Limited
Level 3, 678 Victoria Street
Richmond, VIC, 3121

T: +61 1800 337 041 | F: +61 1800 803 578

North America Region

Ansell Healthcare Products LLC
111 Wood Avenue South, Suite 210
Iselin, NJ 08830 USA

T: +1 800 800 0444 | F: +1 800 800 0445

Europe, Middle East and Africa (EMEA) Region

Ansell Healthcare Europe NV
Riverside Business Park,
Boulevard International, 55
1070 Brussels, Belgium

T: +32 2 528 74 00 | F: +32 2 528 74 01

Latin America and Caribbean Region

Ansell Commercial México S.A. de C.V.
Bvd. Bernardo Quintana No. 7001-C, Q7001 Torre II.
Suites 1304, 1305 y 1306. Col. Centro Sur, C.P. 76079
Querétaro, Qro. México

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